

Keypad Overview for:

ML /MP 155, 285, 305C, 405 MF 125, 285 (Keypad 2)

Alarm History - keys

Display of information on alarms:

- max or min temperature
- average temperature during alarm
- alarm duration

- "WARM" - alarm
- "COLD" - alarm

Alarm LED

Flashing in temperature alarm situation

Parameter keys

- Access parameter mode (for 5 seconds)
- next parameter
- previous parameter
- Change parameter value
- Confirm parameter change

Power fail
Flashing in Power fail alarm situation



Main switch

O = off
I = on

- Power OK

Reset-key / Mute-key

temporarily switches off buzzer in alarm function

Light key

- LED-Test, when pressed during 5 seconds
- Temperature display during power-fail